

# Manual

innuPREP DNase I

## Content

<b>1</b>	<b>Product and order number .....</b>	<b>I</b>
<b>2</b>	<b>Storage conditions .....</b>	<b>I</b>
<b>3</b>	<b>Description.....</b>	<b>II</b>
3.1	Unit definition .....	II
3.2	Quality control .....	II
<b>4</b>	<b>Delivered components.....</b>	<b>III</b>
4.1	innuPREP DNase I .....	III
4.2	buffers .....	III
<b>5</b>	<b>Protocol.....</b>	<b>IV</b>
5.1	Removal of DNA from RNA preparations.....	IV
5.2	Ethanol precipitation .....	V
5.3	Hints and notes .....	V
<b>6</b>	<b>Related products .....</b>	<b>VI</b>

## Section I

---

### 1 Product and order number

<b>Name</b>	<b>Amount</b>	<b>Order-no.</b>
innuPREP DNase I	5.000 Units	845-KS-5210005
innuPREP DNase I	10.000 Units	845-KS-5210010

### 2 Storage conditions

Store innuPREP DNase I at -20 °C in a freezer with constant temperature conditions.

If stored as recommended the innuPREP DNase I will remain stable at least for 12 months.

### **3 Description**

The innuPREP DNase I has been developed for efficient digestion of DNA. For use whenever the presence of small quantities of contaminating DNA may affect the performance of assays, including RT-PCR – innuPREP DNase I demonstrates high nuclease and is certified free from RNase activity.

innuPREP DNase I is an endonuclease which causes breaks of single- and double-stranded DNA resulting mono and oligo nucleotides. DNase I has no influence to RNA which will remain intact. The enzyme is used to remove unmeant co-purified genomic DNA from RNA preparations.

innuPREP DNase I acts independently of DNA sequence to cleave single-stranded, double-stranded DNA and chromatin to produce 5'-phosphate terminated polynucleotides with a 3' positioned free hydroxyl group.

#### **3.1 Unit definition**

One Kunitz unit is the amount of enzyme which is needed for increase of absorption (A<sub>260 nm</sub>) by 0.001 per minute and milliliter at 25 °C and pH 5.0.

#### **3.2 Quality control**

No RNase activity. No Inhibitors.

Functionality tested by digesting template DNA.

## 4 Delivered components

### 4.1 innuPREP DNase I

**Concentration:** 1 Kunitz Units/ $\mu$ l

**DNase I in storage buffer** 

The enzyme is supplied in:

50 mM Tris-HCl (pH 7,5), 10 mM  $\text{CaCl}_2$  and 50% glycerol.

### 4.2 buffers

The innuPREP DNase I (1 KU/ $\mu$ l) is provided with the following buffers:

**DNase I Reaction Buffer (10x)** 

100 mM Tris-HCl (pH 7.5 at 25 °C), 25 mM  $\text{MgCl}_2$  and 1 mM  $\text{CaCl}_2$ .

**EDTA Solution** 

50 mM EDTA.

## 5 Protocol

- Recommended protocols for digestion in combination with the:
  - innuPREP RNA Mini Kit
  - innuPREP RNA MIDI Direct Kit
  - innuPREP Blood RNA Kit
  - innuPREP Blood RNA MIDI Direct Kit
  - innuPREP Plant RNA Kit
  - innuPREP Micro RNA Kit
  - innuSPEED Tissue RNA Kit
  - innuSPEED Plant RNA Kit
  - innuPREP total RNA Kit
  
- After purification of the nucleic acid it could necessary to perform a DNase I step with the eluates of nucleic acids.

### 5.1 Removal of DNA from RNA preparations

- Prepare the following reaction mix in a DNase-/RNase-free reaction tube:

---

1  $\mu$ l innuPREP DNase I (1 Kunitz Units/ $\mu$ l) / 1  $\mu$ g RNA

---

x  $\mu$ l DNase I Reaction Buffer (10x)

---

add with RNase-free water to the final reaction volume

- Incubate the reaction mix for 30 min at 37 °C.
- To inactivate the enzyme add 1  $\mu$ l 50 mM EDTA or EGTA per 10  $\mu$ l reaction mix and incubate the mixture for 10 min at 65 °C.

## 5.2 Ethanol precipitation

- Depending on the downstream application an ethanol precipitation can be reasonable:
- Therefore add 2.5 x volume 100 % Ethanol (pre-cooled at -20 °C) to the RNA solution, mix thoroughly and incubate the mix for minimum 10 min at -20 °C.
- Centrifuge the mixture and remove the supernatant and wash the RNA pellet once with 70 % ethanol by vortexing and subsequent centrifugation for 8 minutes at 7.500 x g (4 °C).
- At the end of the procedure briefly air-dry the pellet for 5 minutes. Dissolve the RNA pellet in DEPCdH<sub>2</sub>O or RNase-free buffer by vortexing or by passing it a few times through a pipette tip.

## 5.3 Hints and notes

- Don't vortex DNase I, mix only by inverting of the tube carefully.
- Use only freshly prepared reaction mix.

## Section VI

### 6 Related products

Product	Amount	Order Number
innuPREP RNA Mini Kit	50 rxn	845-KS-2040050
	250 rxn	845-KS-2040250
innuPREP Blood RNA Kit	50 rxn	845-KS-2010050
	250 rxn	845-KS-2010250
innuPREP Plant RNA Kit	50 rxn	845-KS-2060050
	250 rxn	845-KS-2060250
innuPREP Micro RNA Kit	50 rxn	845-KS-2030050
	250 rxn	845-KS-2030250
innuPREP FFPE total RNA Kit	50 rxn	845-KS-2050050
	250 rxn	845-KS-2050250
innuSPEED Tissue RNA Kit	50 rxn	845-KS-2540050
	250 rxn	845-KS-2540250
innuSPEED Plant RNA Kit	50 rxn	845-KS-2560050
	250 rxn	845-KS-2560250

Publication No.: HB\_KS-5210\_e\_151012

This documentation describes the state at the time of publishing. It needs not necessarily agree with future versions. Subject to change!

*Expression and further use permitted with indication of source. © 2015 Analytik Jena AG, AJ Innuscreen GmbH*

**Manufacturer:**

AJ Innuscreen GmbH

Robert-Rössle-Straße 10  
13125 Berlin

**Distribution/Publisher:**

Analytik Jena AG

Phone +49 (0) 36 41 / 77-94 00  
Fax +49 (0) 36 41 / 77-76 77 76

Konrad-Zuse-Straße 1  
07745 Jena/ Germany  
www.bio.analytik-jena.com  
lifescience@analytik-jena.com

