

Instructions for Use

Life Science Kits & Assays

innuPREP Virus DNA/RNA Kit – KFml

Order No.:

845-KF-4615015 15 reactions

845-KF-4615250 250 reactions

845-KF-4615750 750 reactions

Publication No.: HB_KF-4615_e_180212

This documentation describes the state at the time of publishing.
It needs not necessarily agree with future versions. Subject to change!

Print-out and further use permitted with indication of source.

© Copyright 2018, Analytik Jena AG, AJ Innuscreen GmbH

Manufacturer:

AJ Innuscreen GmbH
Robert-Rössle-Straße 10
13125 Berlin
Made in Germany!

Distribution/Publisher:

Analytik Jena AG
Konrad-Zuse-Straße 1
07745 Jena · Germany

Phone +49 3641 77 9400
Fax +49 3641 77 767776
www.analytik-jena.com
info@analytik-jena.com

Contents

1	Introduction.....	3
1.1	Intended use.....	3
1.2	Notes on the use of this manual	4
2	Safety precautions	5
3	General notes and safety recommendations on handling RNA.....	6
4	Storage conditions	7
5	Functional testing and technical assistance	8
6	Product use and warranty	8
7	Kit components	9
8	Recommended steps before starting	11
9	Carrier Mix.....	11
9.1	Storage conditions and handling	11
9.2	Preparation of Lysis Solution RL/Carrier Mix.....	12
10	GHS Classification	13
10.1	Hazard phrases	13
10.2	Precaution phrases	14
10.3	EU hazard statements	14
11	Product specifications.....	15
12	Lysis protocols for isolation of viral DNA/RNA from different starting materials.....	16
12.1	Protocol 1: Isolation of viral DNA and RNA from 200 µl cell-free body fluids (serum, plasma, cerebrospinal fluid, liquor) ..	16
12.2	Protocol 2: Isolation of viral DNA/RNA from cell culture supernatants	16
12.3	Protocol 3: Isolation of viral DNA/RNA from swab samples ...	17
12.4	Protocol 4: Isolation of viral DNA/RNA from tissue biopsies ..	18
12.5	Protocol 5: Isolation of viral DNA/RNA stool samples (tested for Norovirus extraction)	19
13	Preliminary steps of the KingFisher ml.....	21

13.1	Filling of KingFisher Tube Strips	21
13.2	Loading the sample	22
14	Automatic processing of the KingFisher ml.....	22
15	Troubleshooting.....	23
16	Related Products	24

1 Introduction

1.1 Intended use

The innuPREP Virus DNA/RNA Kit - KFml has been designed for isolation of viral DNA/RNA from different kinds of starting material. The kit contains a Carrier Mix with Carrier RNA as well as an internal extraction controls for DNA and RNA. The Internal Control DNA or RNA can be detected by real-time PCR using the corresponding assays.

The extraction procedure is based on a new kind of chemistry (patent pending). The procedure combines lysis of starting material with subsequent binding of viral DNA/RNA on surface modified magnetic particles. After washing steps the viral DNA/RNA is eluted from the magnetic particles by using water.

The extraction procedure takes place on the magnetic particle processor KingFisher ml. Extraction chemistry and extraction protocol are optimized to get maximum yield of RNA.



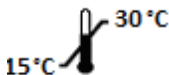





CONSULT INSTRUCTION FOR USE



This package insert must be read carefully prior to use. Package insert instructions must be followed accordingly. Reliability of results cannot be guaranteed if there are any deviations from the instructions in this package insert.

1.2 Notes on the use of this manual

For easy reference and orientation, the manual uses the following warning and information symbols as well as the shown methodology:

Symbol	Information
	REF Catalogue number.
	Content Contains sufficient reagents for <N> tests.
	Storage conditions Store at room temperature, unless otherwise specified.
	Consult instructions for use This information must be observed to avoid improper use of the kit and the kit components.
	Expiry date
	Lot number The number of the kit charge.
	Manufactured by Contact information of manufacturer.
	For single use only Do not use components for a second time.
	Note / Attention Observe the notes marked in this way to ensure correct function of the device and to avoid operating errors for obtaining correct results.

2 Safety precautions

NOTE

Read through this chapter carefully prior to use to guarantee your own safety and a trouble-free operation.

Follow all the safety instructions explained in the manual, as well as all messages and information, which are shown.

All due care and attention should be exercised in handling the materials and reagents contained in the kit. Always wear gloves while handling these reagents and avoid any skin contact! In case of contact, flush eyes or skin with a large amount of water immediately.



FOR SINGLE USE ONLY!

This kit is made for single use only!

ATTENTION!

Don't eat or drink components of the kit!

The kit is designed to be handled by educated personnel in a laboratory environment!

If the buffer bottles are damaged or leaking, wear gloves and protective goggles when discarding the bottles in order to avoid any injuries. This kit is to be used with potential infectious human samples. Therefore, all liquid waste must be considered as potentially infectious and must be handled and discarded according to local safety regulation.

Please observe the federal, state and local safety and environmental regulations. Follow the usual precautions for applications using extracted nucleic acids. All materials and reagents used for DNA or RNA isolation should be free of DNases or RNases.

ATTENTION!

Do not add bleach or acidic components to the waste after sample preparation!

NOTE

Emergency medical information in English and German can be obtained 24 hours a day from:

Poison Information Center, Freiburg / Germany

Phone: +49 (0)761 19 240.

For more information, please ask for the material safety data sheets (MSDS's).

3 General notes and safety recommendations on handling RNA

RNA is far less stable than DNA. It is very sensitive to degradation by endogenous RNases in the biological material and exogenous RNases which are permanently present everywhere in the lab. To achieve satisfactory qualitative and quantitative results in RNA preparations, contaminations with exogenous RNases have to be reduced to a minimum in accordance with the following recommendations:

- Always wear latex or vinyl gloves while handling reagents and RNA samples to prevent RNase contaminations from surface of the skin or from dusty laboratory equipment.
- Change gloves frequently and keep tubes closed.
- Keep isolated RNA on ice.
- Reduce preparation time as much as possible.
- Use only sterile, disposable polypropylene tubes throughout the procedure (these tubes are generally RNase-free.)
- Non-disposable plastic ware should be treated before use to ensure that it is RNase-free. Plastic ware should be thoroughly rinsed with 0.1 M NaOH, 1 mM EDTA followed by RNase-free water. You can also take chloroform-resistant plastic ware rinsed with chloroform to inactivate RNases.

- All glassware should be treated before use to ensure that it is RNase-free. Glassware should be cleaned with detergent, thoroughly rinsed and oven baked at 240 °C for four hours or more before use. Autoclaving will not inactivate RNase activity completely. Oven baking inactivates RNases and ensures that no other nucleic acids (such as plasmid DNA) are present on the surface of the glassware. You can also clean glassware with 0.1 % DEPC (diethyl pyrocarbonate). The glassware has to be immersed in 0.1 % DEPC solution for 12 hours at 37 °C followed by autoclaving or heating to 100 °C for 15 minutes to remove residual DEPC.
- Electrophoresis tanks should be cleaned with detergent solution (e.g. 0.5 % SDS), thoroughly rinsed with RNase-free water, rinsed with ethanol and finally allowed to dry.
- All buffers have to be prepared with DEPC-treated RNase-free water.
- Avoid handling bacterial cultures, cell cultures or other biological sources of RNases in the same lab where the RNA purification will be performed.
- Do not use equipment, glassware and plastic ware employed for other applications which might introduce RNase contaminations in the RNA isolation.

4 Storage conditions

Store lyophilized **Carrier Mix** at -22 °C to -18 °C.

All other components of the innuPREP Virus DNA/RNA Kit - KFml should be stored dry at room temperature (15 °C to 30 °C). When stored at room temperature, the kit is stable until the expiration date printed on the label on the kit box.

Before every use make sure that all components have room temperature. If there are any precipitates within the provided solutions dissolve these precipitates by careful warming.

For further information see chapter "Kit components" p. 9.

5 Functional testing and technical assistance

The Analytik Jena AG guarantees the correct function of the kit for applications as described in the manual. This product has been produced and tested in an ISO 13485 certified facility.

We reserve the right to change or modify our products to enhance their performance and design. If you have any questions or problems regarding any aspects of the innuPREP Virus DNA/RNA Kit - KFml or other Analytik Jena AG products, please do not hesitate to contact us. For technical support or further information in Germany please dial +49 36 41 / 77 94 00. For other countries please contact your local distributor.

6 Product use and warranty

The kit is not designed for the usage of other starting materials or other amounts of starting materials than those, referred to in the manual (→ "Intended use" p. 3) (→ "Product specifications" p. 15). Since the performance characteristics of Analytik Jena AG kits have just been validated for the application described above, the user is responsible for the validation of the performance of Analytik Jena AG kits using other protocols than those described below. Analytik Jena AG kits may be used in clinical diagnostic laboratory systems after the laboratory has validated the complete diagnostic system as required by CLIA' 88 regulations in the U.S. or equivalent regulations required in other countries.

All products sold by the Analytik Jena AG are subjected to extensive quality control procedures and are warranted to perform as described when used correctly. Any problems should be reported immediately.

NOTE

The kit is for research use only !

7 Kit components




IMPORTANT

Store lyophilized Carrier Mix at -22 °C to -18 °C.



STORAGE CONDITIONS

All other components are stored at room temperature.

	 15	 250	 750
REF	845-KF-4615015	845-KF-4615250	845-KF-4615750
MAG Suspension	1 ml	2 x 9 ml	5x 9 ml
Carrier Mix	For 1 x 1.25 ml working solution	For 4 x 1.25 ml working solution	For 10 x 1.25 ml working solution
RNase-free Water	2 ml	4x 2 ml	10x 2 ml
Lysis Solution RL	10 ml	140 ml	2x 200 ml
Binding Solution RBS	10 ml	125 ml	2x 180 ml
Washing Solution HS (conc.)	5 ml	70 ml	3x 70 ml
Washing Solution LS (conc.)	6 ml	2x 50 ml	2x 130 ml
RNase-free Water	2 ml	2x 25 ml	4x 25 ml
KingFisher Tip Combs	3	50	150
KingFisher Tube Strips	15	250	750
Elution Tubes	15	5x 50	15x 50
Manual	1	1	1

Kit components

Initial steps

Carrier Mix

Dissolve Carrier Mix by addition of 1.25 ml RNase-free Water, mix thoroughly by pipetting up and down!

Washing Solution HS

Add 5 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

Washing Solution LS

Add 24 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

Carrier Mix

Dissolve Carrier Mix by addition of 1.25 ml RNase-free Water, mix thoroughly by pipetting up and down!

Washing Solution HS

Add 70 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

Washing Solution LS

Add 200 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

Carrier Mix

Dissolve Carrier Mix by addition of 1.25 ml RNase-free Water, mix thoroughly by pipetting up and down!

Washing Solution HS

Add 70 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

Washing Solution LS

Add 560 ml of 96-99.8 % ethanol to the bottle and mix thoroughly. Keep the bottle always firmly closed!

COMPONENTS NOT INCLUDED IN THE KIT

- 96–99.8 % ethanol, non-denatured or methylated 1.5 ml tubes
- 1 x PBS (137 mM NaCl, 2.7 mM KCl, 10 mM Na₂HPO₄, 1.8 mM KH₂PO₄)
- Physiological saline

8 Recommended steps before starting

- Ensure that the **Washing Solution HS** and **Washing Solution LS** have been prepared according to the instruction (→ "Kit components" p. 9).
- Ensure that the **Carrier Mix** and **Lysis Solution RL / Carrier Mix** have been prepared according to the instruction (→ "Carrier Mix" p. 11).
- Centrifugation steps should be carried out at room temperature.
- Avoid freezing and thawing of starting material.

9 Carrier Mix

9.1 Storage conditions and handling

The Carrier Mix contains a carrier RNA and an internal control RNA and DNA (IC RNA and IC DNA).

- Add dissolved Carrier Mix to Lysis Solution RL immediately.
- Unused Carrier Mix should be kept frozen at -18 °C to -22 °C.
- Do not freeze and thaw the Carrier Mix more than 3 times.
- Mixture of Lysis Solution RL and Carrier Mix is stable for 7 days at 4 °C.
- Internal control RNA or DNA can be detected by Real-time PCR using the corresponding assays, as shown below:

Name	Amount	Order-no.
Detection of IC RNA		
innuDETECT Internal Control RNA Assay	100 rxn	845-ID-0006100
Detection of IC DNA		
innuDETECT Internal Control DNA Assay	100 rxn	845-ID-0007100
Detection of IC RNA and IC DNA		
innuDETECT Internal Control DNA/RNA Assay	100 rxn	845-ID-0008100

9.2 Preparation of Lysis Solution RL/Carrier Mix

1. Add 1.25 ml RNase-free Water to each tube **Carrier Mix**.
2. Mix thoroughly by pipetting up and down!
3. After the preparation of **Carrier Mix** stock solution prepare the mixture of **Lysis Solution RL/Carrier Mix** as described in the following table:

Component	1 samples	5 samples	15 samples	n samples
Lysis Solution RL	0.5 ml	2.5 ml	7.5 ml	500 µl x n samples
Carrier Mix	16.5 µl	82.5 µl	247.5 µl	16.5 µl x n samples

NOTE

If customized Extraction Controls should be used, please add these components to the **Lysis Solution RL/Carrier Mix**!

10 GHS Classification

Component	Hazard contents	GHS Symbol	Hazard phrases	Precaution phrases	EUH
Lysis Solution RL	Guanidinium thiocyanate 25-50 %	The GHS symbols for this component are Corrosive (C) and Hazardous (Xn), both shown in red diamond shapes.	302, 314, 412	101, 102, 103, 260,303+361+353, 305+351+338, 310, 405, 501	032
Danger					
Binding Solution RBS	Guanidinium thiocyanate 25-50 %	The GHS symbols for this component are Corrosive (C) and Hazardous (Xn), both shown in red diamond shapes.	302, 314, 412	101, 102, 103, 260,303+361+353, 305+351+338, 310, 405, 501	032
Danger					
Washing Solution HS (conc.)	Guanidinium thiocyanate 50-100 %	The GHS symbols for this component are Hazardous (Xn) and Corrosive (C), both shown in red diamond shapes.	302, 314, 412	101, 102, 103, 260,303+361+353, 305+351+338, 310, 405, 501	032
Danger					

10.1 Hazard phrases

- 302 Harmful if swallowed.
- 314 Causes severe skin burns and eye damage.
- 412 Harmful to aquatic life with long lasting effects.

10.2 Precaution phrases

- 101 If medical advice is needed, have product container or label at hand.
- 102 Keep out of reach of children.
- 103 Read label before use.
- 260 Do not breathe dust/fume/gas/mist/vapors/spray.
- 310 Immediately call a POISON CENTER/doctor.
- 405 Store locked up.
- 501 Dispose of contents/container in accordance with local/regional/national/international regulations.
- 303+361+353 IF ON SKIN (or hair): Take off immediately all contaminated clothing. Rinse skin with water/shower.
- 305+351+338 IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

10.3 EU hazard statements

- 032 Contact with acids liberates very toxic gas.

11 Product specifications

1. Starting material:

- Cell-free body fluids (serum, plasma, cerebrospinal fluid, liquor); max. 200 µl sample volume
- Cell culture supernatants; max. 200 µl sample volume
- Swab samples ; max. 200 µl sample volume
- Stool samples ; max. 200 µl sample volume
- Tissue biopsies (1–5 mg)

12 Lysis protocols for isolation of viral DNA/RNA from different starting materials

12.1 Protocol 1: Isolation of viral DNA and RNA from 200 µl cell-free body fluids (serum, plasma, cerebrospinal fluid, liquor)

1. Transfer 300 µl Lysis Solution RL/Carrier Mix into a 1.5 ml Tube and add 200 µl of the sample.
2. Close the cap and vortex the 1.5 ml Tube for 10 seconds.
3. Place the 1.5 ml Tube into a thermal mixer and incubate under continuous shaking for 15 minutes at room temperature.

NOTE

Alternative the 1.5 ml Tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds).

12.2 Protocol 2: Isolation of viral DNA/RNA from cell culture supernatants

1. Transfer 300 µl Lysis Solution RL/Carrier Mix into a 1.5 ml Tube and add 200 µl of the cell culture supernatant (cell culture medium).
2. Close the cap and vortex the 1.5 ml Tube for 10 seconds.
3. Place the 1.5 ml Tube into a thermal mixer and incubate under continuous shaking for 15 minutes at room temperature.

NOTE

Alternative the 1.5 ml Tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds).

12.3 Protocol 3: Isolation of viral DNA/RNA from swab samples

1. Place the swab into a 1.5 ml reaction tube and add **500 µl Lysis Solution RL/Carrier Mix**. Vortex shortly!

NOTE

To get maximum yield of viral nucleic acids it is essential to leave the swab in the reaction tube during the complete lysis time. It is possible to cut the shaft of the swab, so that you can close the cap of the reaction tube. It is also possible to perform the lysis step with opened cap. The removal of the swab from the reaction tube ahead of time will lead to a dramatically reduced final yield!

2. Place the reaction tube into a thermal mixer and incubate under continuous shaking for 15 minutes at room temperature.

NOTE

Alternative the 1.5 ml Tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds).

3. After lysis time carefully squeeze out the swab on the wall of the tube and discard the swab.

12.4 Protocol 4: Isolation of viral DNA/RNA from tissue biopsies

1. Transfer 500 µl Lysis Solution RL/Carrier Mix into a 1.5 ml reaction tube and add about 1–5 mg of the tissue biopsy
2. Close the cap and vortex the reaction tube for 10 seconds.
3. Place the reaction tube into a thermal mixer and incubate under continuous shaking for 30 minutes at room temperature. Lysis time can be increased up to lysis of starting material is complete (60 minutes).

NOTE

Alternative the reaction tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds). A longer lysis time can lead to a reduced yield and quality of some viral DNA/RNA.

4. After lysis centrifuge the reaction tube at max. speed for 1 minute to spin down unlysed material and follow the manual exactly for the next steps.

12.5 Protocol 5: Isolation of viral DNA/RNA stool samples (tested for Norovirus extraction)

Alternative 1: Solid stool sample

1. Transfer about **0.05 – 0.1 g** of the **stool sample** into a 1.5 ml reaction tube and add **200 µl PBS** (not included in the kit).
2. Vortex the sample for 5 seconds and centrifuge it at max. speed for 3 minutes.
3. Transfer respectively **300 µl Lysis Solution RL/Carrier Mix** into an reaction tube and add the **clarified supernatant** of the stool sample from step 2.
4. Close the cap and vortex the reaction tube for 10 seconds.
5. Place the reaction tube into a thermal mixer and incubate under continuous shaking for 15 minutes at room temperature.

NOTE

Alternative the 1.5 ml Tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds).

Alternative 2: Homogenized stool sample

1. Centrifuge the homogenized stool sample to spin down still remaining particles.
2. Transfer 300 µl Lysis Solution RL/Carrier Mix into a reaction tube and add the 200 µl of the clarified supernatant of the stool sample from step 1.
3. Close the cap and vortex the reaction tube for 10 seconds.
4. Place the reaction tube into a thermal mixer and incubate under continuous shaking for 15 minutes at room temperature.

NOTE

Alternative the 1.5 ml Tube can be mixed by vortexing during the lysis (each 5 minutes for 5 seconds).

13 Preliminary steps of the KingFisher ml

13.1 Filling of KingFisher Tube Strips

NOTE

During sample lysis pre-fill the tubes of the KingFisher Tube Strips with the following buffers respectively.

IMPORTANT

It is important to mix the **MAG Suspension** by vigorous shaking or vortexing before use (approx. 30 seconds)!

Filling of KingFisher Tube Strips

Tube A	450 µl	Binding Solution RBS
	50 µl	MAG Suspension
Tube B	500 µl	Washing Solution HS
Tube C	800 µl	Washing Solution LS
Tube D	800 µl	Washing Solution LS
Tube E	120 µl	RNase-free Water

13.2 Loading the sample

NOTE

The following step will be done after the sample lysis!

After lysis (in case of the biopsy sample after centrifugation) transfer the lysed sample (approx. 500 µl) into the **Tube A** of the KingFisher Tube Strip.

The final volume has to be 1 ml (lysed sample and **Binding Solution RBS** and **MAG Suspension**).

14 Automatic processing of the KingFisher ml

1. Place the filled KingFisher Tube Strips into the KingFisher system on the right position!
2. Place the KingFisher Tip Combs onto the magnetic track!
3. Start the program "INNU_ViralRNA_KFml"!

NOTE

After finishing the extraction protocol, the **Tube E** of the Tube Strip contains the extracted RNA. Store the RNA under adequate conditions. We recommend to store the extracted RNA at $-80\text{ }^{\circ}\text{C}$.

15 Troubleshooting

Problem / probable cause	Comments and suggestions
Low amount of extracted viral RNA/DNA	
Insufficient lysis of starting material.	<p>Ensure to use the correct mixture of Lysis Solution RL and Carrier Mix.</p> <p>Ensure to incubate at room temperature for sample lysis.</p>
Inadequate extraction.	<p>Inhibiting substances in starting material. Please use the kit only for samples that match the requirements declared in "Product specifications".</p> <p>Use Internal Controls for verification of extraction procedure.</p>

16 Related Products

Name	Amount	Order No.
Detection of Internal Controls		
Detection of IC DNA		
innuDETECT Internal Control DNA Assay	100 rxn	845-ID-0006100
Detection of IC RNA		
innuDETECT Internal Control RNA Assay	100 rxn	845-ID-0007100
Detection of IC DNA and IC RNA		
innuDETECT Internal Control DNA/RNA Assay	100 rxn	845-ID-0008100
Products for PCR & Electrophoresis		
innuTaq DNA Polymerase (5 U/μl)	500 U	845-EZ-1000500
innuTaq Hot-A DNA Polymerase (5 U/μl)	500 U	845-EZ-3000500
innuMIX Standard PCR MasterMix	100 rxn	845-AS-1700100
	200 rxn	845-AS-1700200
innuMIX Green PCR MasterMix	100 rxn	845-AS-1400100
	200 rxn	845-AS-1400200

Headquarters

Analytik Jena AG
Konrad-Zuse-Str. 1
07745 Jena · Germany

Phone +49 3641 77 70
Fax +49 3641 77 9279
info@analytik-jena.com
www.analytik-jena.com

Pictures: Analytik Jena AG
Subject to changes in design and scope of delivery as well as further technical development!