
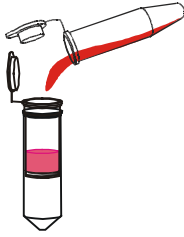
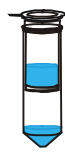




innuPREP DNA Micro Kit

Protocol 1: DNA isolation from small tissue samples or biopsies

- Recommended steps before starting
- Heat thermal mixer or water bath (50 °C)
 - Prepare Washing Solution HS, Washing Solution MS and Proteinase K according to the instruction

1. Starting material		<ul style="list-style-type: none"> ▪ Max. 5 mg
2. Lysis		<ul style="list-style-type: none"> ▪ Add 200 µl Lysis Solution TLS <u>and</u> 20 µl Proteinase K ▪ Vortex: 5 sec ▪ Incubation: 50 °C, 30 min ▪ 10.000 x g (~12.000 rpm): 1 min ▪ Transfer supernatant; 1.5 ml tube
3. Optional: RNA removal		<ul style="list-style-type: none"> ▪ 4 µl 100 mg/ml RNase A; vortex ▪ Incubation: 5 min @ RT
4. Binding of DNA		<ul style="list-style-type: none"> ▪ Add 200 µl Binding Solution TBS ▪ Vortex: 15 sec ▪ Add Spin Filter to Receiver Tube ▪ Add sample to Spin Filter ▪ 10.000 x g (~12.000 rpm): 1 min
5. Washing		<ul style="list-style-type: none"> ▪ Add 400 µl Washing Solution HS ▪ 10.000 x g (~12.000 rpm): 30 sec ▪ Add 750 µl Washing Solution MS ▪ 10.000 x g (~12.000 rpm): 30 sec
6. Remove Ethanol		<ul style="list-style-type: none"> ▪ Discard filtrate ▪ Add Spin Filter to Receiver Tube ▪ Centrifuge: max speed, 2 min
7. Elution		<ul style="list-style-type: none"> ▪ Add Spin Filter to an Elution Tube ▪ Add 50 – 100 µl Elution Buffer ▪ Incubation: 1 min @ RT ▪ 6.000 x g (~8.000 rpm): 1 min

Order No.: 845-KS-1011010 10 reactions
845-KS-1011050 50 reactions

This documentation describes the state at the time of publishing. It needs not necessarily agree with future versions. Subject to change!

Expression and further use permitted with indication of source. © 2018 Analytik Jena AG, AJ Innuscreen GmbH

Manufacturer:

AJ Innuscreen GmbH

Robert-Rössle-Straße 10
13125 Berlin

Distribution/Publisher:

Analytik Jena AG

Konrad-Zuse-Straße 1
07745 Jena · Germany
www.analytik-jena.com
info@analytik-jena.com

Phone +49 36 41 77 9400
Fax +49 36 41 77 767776